

**ROWLEY**  
**BIOCHEMICAL**  
**INC.**

DANVERS INDUSTRIAL PARK  
10 ELECTRONICS AVENUE,  
DANVERS MA 01923  
TEL: 978-739-4883 FAX: 978-739-5640  
www.rowleybio.com

---

**L-759**      **GILL'S HEMATOXYLIN #1**  
**FOR CYTOLOGY**

**FIXATION:** Wet-fix in 95% Ethyl alcohol.

**STAINING PROCEDURE:**

1. Ethyl alcohol, 95% for 15 minutes.
2. Two changes distilled water, 10 dips each.
3. **Gill's Hematoxylin # 1** (L-759-1) for 2 minutes.
4. Two changes distilled water, 10 dips each.
5. **Scott's Tap Water Substitute** (L-759-4) for 1 minute.
6. Two changes distilled water, 10 dips each.
7. Ethyl alcohol, 95% 10 dips and counterstain.

**RESULTS:**

Chromatin              Blue to Blue Black

**REFERENCES:**

Gill, G.W., Acta Cytol., 18:300-311, 1974