

**L-759 GILL'S HEMATOXYLIN, # 2 & # 3,
FOR HISTOLOGY**

Fixation: Fixative of choice.

Sections: Paraffin

Staining:

1. Deparaffinize and hydrate to distilled water.
2. Stain in **Gill's Hematoxylin # 2** (L-759-2) for 3 minutes. If a more intense hematoxylin stain is desired stain in **Gill's Hematoxylin # 3** (L-759-3),
3. Wash in two changes of distilled water, 10 dips each.
4. Blue sections in **Scott's Tap Water Substitute** (L-759-4) for 1 minute.
5. Wash thoroughly and counterstain.

RESULTS:

Chromatin	Blue to Blue Black
-----------	--------------------

REFERENCES:

Gill, G.W., Acta Cytol., 18:300-311, 1974.