

**ROWLEY
BIOCHEMICAL
INC.**

**DANVERS INDUSTRIAL PARK
10 ELECTRONICS AVENUE,
DANVERS MA 01923
TEL: 978-739-4883 FAX: 978-739-5640**

L-760, CONGO RED

STAINING:

1. Well washed sections or slides stained with an Alum Hematoxylin.
Staining the sections or slides in **Congo Red Staining Solution** (L-760-1), for 1 minute.
2. Rinse in water.
3. Dehydrate , clear and mount.

Stain Results:

Nuclei	Blue
Cytoplasm	Pink
Collagen	Reticulocytes

References:

Clark, G.: Staining Procedures, Williams & Wilkins Co., Baltimore,
3rd Ed., c. 1973, p. 37.

Conn, H.J.: Biological Stains, 6th ed., 1953 Williams and Wilkins Co., Baltimore.